



KAPA SYBR® FAST One-Step qRT-PCR Kits

KAPA SYBR® FAST One-Step qRT-PCR Kits contain M-MuLV Reverse Transcriptase, RNase Inhibitor and a novel DNA Polymerase engineered via molecular evolution. The kit is optimized for rapid one-step, one-tube RNA quantification.

KAPA SYBR® FAST One-Step qRT-PCR Kits are designed to perform optimally in stringent qRT-PCR reaction conditions, exhibiting dramatic improvements in sensitivity, specificity, reaction efficiency, and fluorescence.




Kits are ideally suited for:

- Gene expression analysis
- Microarray validation
- Gene knockdown validation
- RNAi and miRNA research

Reduce experimental variation and contamination with a convenient one-step qRT-PCR protocol.

Superior signal and reaction efficiency


The improved speed, processivity and robustness of KAPA SYBR® FAST One-Step qRT-PCR kits results in consistently high amplification efficiencies required for accurate relative quantification. To demonstrate the high performance of KAPA SYBR® FAST One-Step qRT-PCR Kit for gene expression analysis, the kit was compared to leading competitor kits across a wide range of input RNA concentrations. The KAPA SYBR® FAST One-Step qRT-PCR Kit achieved higher fluorescence, earlier Ct values and improved amplification efficiency when compared to competitor one-step qRT-PCR kits.

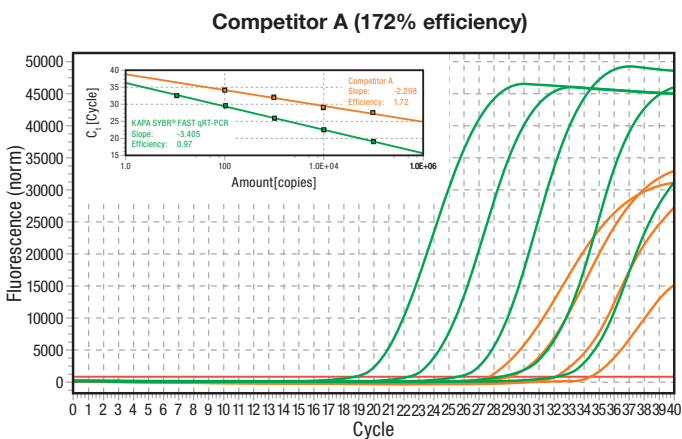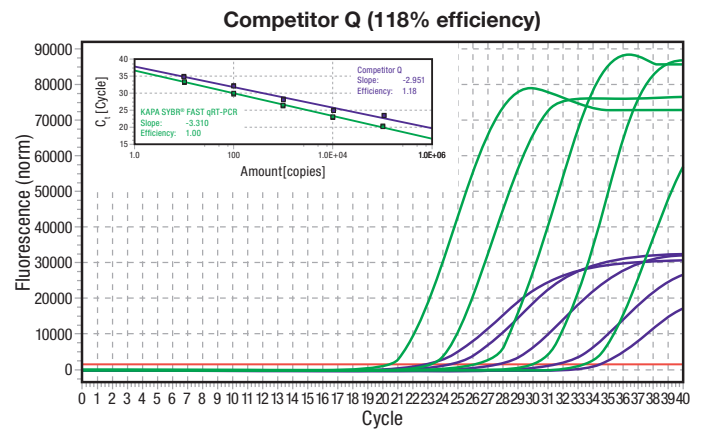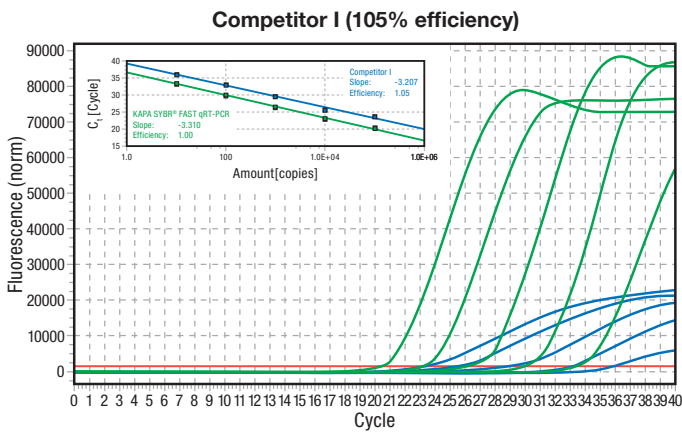

Higher fluorescence, earlier Ct values and improved reaction efficiency. The RRMI gene (94 bp, 45.7% GC) was amplified from log-fold serial dilutions of RNA (100 ng- 10 pg/20 µl reaction), isolated from human placenta, using the KAPA SYBR® FAST One-Step qRT-PCR Kit (green) or Competitor I (blue), Competitor Q (purple), and Competitor A (orange). Linear amplification plots demonstrate earlier Ct values and greater baseline subtracted fluorescence for the RRMI gene with the KAPA SYBR® FAST One-Step qRT-PCR Kit. Calculated reaction efficiencies confirmed consistently high performance is achievable with the KAPA SYBR® FAST One-Step qRT-PCR Kit (RRMI = 100% and 97%). Efficiencies obtained with Competitor I (105%), Competitor Q (118%) and Competitor A (176%) were sub-optimal. Reactions were performed using the competitors recommend protocols on the Eppendorf® MasterCycler™ ep realplex⁴ S real-time PCR cycler.


KAPA SYBR® FAST One-Step qRT-PCR Kits

Robust amplification across a broad range of target lengths

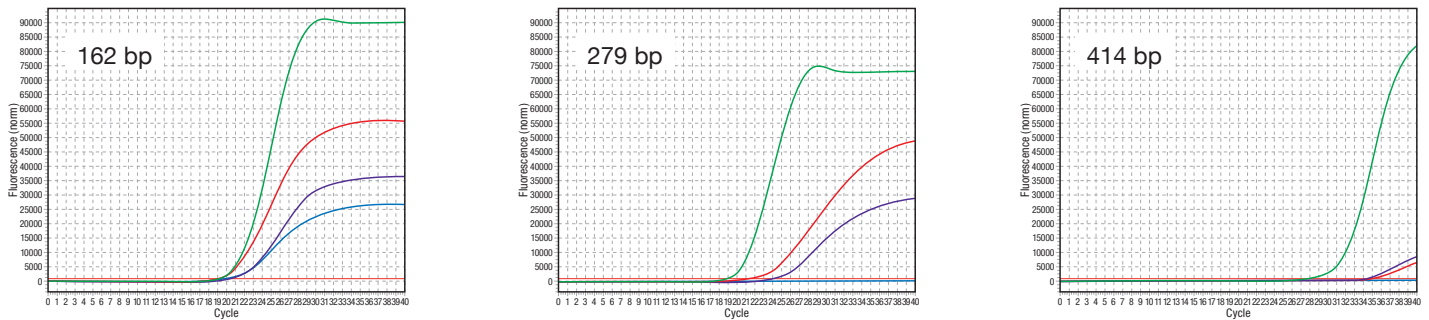

To demonstrate the robust performance of the KAPA SYBR® FAST One-Step qRT-PCR Kit, MS2 bacteriophage RNA targets of increasing length were amplified at a concentration of 0.8 pg/20 µl using the KAPA SYBR® FAST One-Step qRT-PCR Kit (green) or Competitor T (red), Competitor I (blue), and Competitor Q (purple). Increased amplicon length led to a significant decrease in competitor performance as indicated by delayed Ct values or complete reaction failure. cDNA synthesis and qPCR amplification were performed using the competitors recommend protocols on the Eppendorf® MasterCycler™ ep *realplex*⁴ S real-time PCR cycler.

High reproducibility and sensitivity across a broad range of concentrations

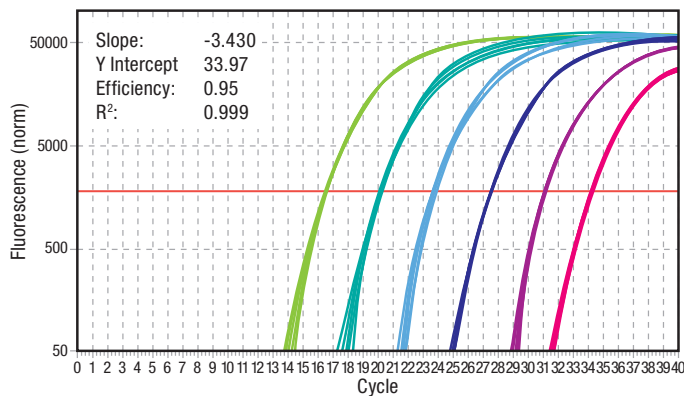

The high reproducibility and sensitivity of the KAPA SYBR® FAST One-Step qRT-PCR Kit enable accurate interrogation of a broad range of RNA template concentrations. A 279 bp fragment was amplified in replicates of six from log-fold serial dilutions of commercial MS2 bacteriophage RNA (80 pg - 0.8 fg/20 µl). cDNA synthesis was performed at 42 °C for 5 min followed by 95 °C for 5 min and 40 cycles of 3 sec denaturation at 95 °C and 30 sec combined annealing/extension at 60 °C. The calculated reaction efficiency of 95% was obtained using all six replicates per data point. Reactions were performed using the competitors recommend protocols on the Eppendorf® MasterCycler™ ep *realplex*⁴ S real-time PCR cycler.



ORDERING INFORMATION

Description	Code	Kit contents
KAPA SYBR® FAST One-Step Universal	KK4650	1 x 1 ml
KAPA SYBR® FAST One-Step Universal	KK4651	1 x 5 ml
KAPA SYBR® FAST One-Step Universal	KK4652	2 x 5 ml
KAPA SYBR® FAST One-Step ABI Prism®	KK4660	1 x 1 ml
KAPA SYBR® FAST One-Step ABI Prism®	KK4661	1 x 5 ml
KAPA SYBR® FAST One-Step ABI Prism®	KK4662	2 x 5 ml
KAPA SYBR® FAST One-Step Bio-Rad iCycler™	KK4670	1 x 1 ml
KAPA SYBR® FAST One-Step Bio-Rad iCycler™	KK4671	1 x 5 ml
KAPA SYBR® FAST One-Step Bio-Rad iCycler™	KK4672	2 x 5 ml
KAPA SYBR® FAST One-Step for Roche LC 480	KK4680	1 x 1 ml
KAPA SYBR® FAST One-Step for Roche LC 480	KK4681	1 x 5 ml
KAPA SYBR® FAST One-Step for Roche LC 480	KK4682	2 x 5 ml



For more information please contact sales@kapabiosystems.com or your local representative.

Purchase of this product includes an immunity from suit under patents specified in the product insert to use only the amount purchased for the purchaser's own internal research. No other patent rights are conveyed expressly, by implication, or by estoppel. Further information on purchasing licenses may be obtained by contacting the Director of Licensing, Applied Biosystems, 850 Lincoln Centre Drive, Foster City, California 94404, USA.

Licensed under U.S. Patent Nos. 5,338,671 and 5,587,287 and corresponding patents in other countries. This product is intended for research use only under patents and patent applications owned by Molecular Probes, Inc. SYBR® is a registered trademark of Molecular Probes, Inc. MasterCycler™, iCycler™, Prism® and LightCycler® are trademarks or registered trademarks of their respective companies.