



## Treatment of Vivaspin® concentrators for improved recovery of low-concentrated protein samples



Application  
Note

#05

#06

#07

#08

#09

### Introduction

With appropriate device size and membrane cut-off selected, Vivaspin® products will typically yield recoveries for the concentrated sample > 90% when the starting sample contains over 0.1 mg/ml protein of interest. Depending on sample characteristics relative to the membrane type used, solute (protein) adsorption on the membrane surface is typically very low (2–10 µg/cm<sup>2</sup>) and in practice not detectable.

This can increase to 20–100 µg/cm<sup>2</sup> when the filtrate is of interest and the sample must pass through the whole internal structure of the membrane. Whilst the relative adsorption to the plastic of the sample container will be proportionately less important than on the membrane, due to the higher total surface area, this can be also be a source of yield loss. Typically, a higher cut-off membrane will bind more than a low molecular weight alternative.

Whenever possible, the smallest MWCO and device size applicable should be chosen. Swinging bucket rotors are preferred to fixed angle rotors. This reduces the surface area of the concentrator that will be exposed to the solution during centrifugation.

An important factor not to be neglected is the thorough recovery of the retentate. Make sure to carefully remove all traces of solution from the sample container and, if feasible, rinse the device after recovering the sample with one or more drops of buffer and then recover again.

The intention of the following “passivation” procedure is to improve recovery of protein samples in the nano- to microgram concentration range by pretreating the device (membrane & plastic). For this purpose a range of solutions are suggested in Table 1.

Type	Concentration
Powdered milk	1% in Arium water
BSA	1% in PBS
Tween 20	5% in Arium water
SDS	5% in Arium water
Triton X-100	5% in Arium water
PEG 3000	5% in Arium water

Table 1: Passivation Solutions

### Passivation procedure for Vivaspin® ultrafiltration concentrators

#### A) Passivation Procedure

1. Wash the concentrators once by filling with Arium water and spin the liquid through according to the respective protocol.
2. Remove residual water thoroughly by pipetting. **Caution: Take care not to damage the membrane with the pipette tip.**
3. Fill concentrators with the blocking solution of choice as given in Table 1.
4. Incubate the filled concentrators at room temperature for at least 2 hours (overnight is also possible except for **Triton X-100 which is not recommended for overnight incubation**).
5. Pour out the blocking solution.
6. Rinse the device 3–4x very thoroughly with Arium water and finally spin through.
7. The “passivated” devices are now ready for use. We recommend comparing different passivation reagents with an untreated device.

## Note

It is necessary to rinse the device thoroughly before each washspin to ensure that traces of passivation compound are removed from the deadstop. Use the device immediately for protein concentration or store it at 4°C filled with Arrium water, to prevent the membrane from drying.

## B) Evaluation of passivation effects (exemplary with BSA)

1. Prepare a 10 µg/ml BSA stock solution e.g. by diluting 90 µl of the 4 mg/ml stock solution in 36 ml 0.1 M sodium borate pH 9.3. Mix well.
2. Fill Vivaspin® 2 devices with 2 ml of this 10 µg/ml BSA solution and close with cap provided.
3. Spin the device in a swing-out rotor at 4,000 × g until the volume is to app. 100 µl.
4. Recover the concentrate and make back up to 2 ml with 0.1 M sodium borate pH 9.3
5. Determine recovered protein concentrations e.g. according to Bradford or BCA assays.


## Results and Discussion

As an example, the effect of milk powder was analysed. It could be shown (Table 2) that the protein recovery of a 10 µg/ml BSA solution could be increased from around 70 to 90%.


If milk powder is not interfering with sample purity and quality, it is a good starting point to improve recovery of diluted sample solutions.

## Protein recovery (10 µg/ml BSA) with Vivaspin® PES 10 kDa after passivation

In another example, detergents were analysed with only 250 and 500 ng BSA (Table 3). BSA recovery declined to 50–30% in untreated devices as the protein concentration was reduced. Significant improvement to 60–90% recovery could be demonstrated when using the passivation strategy. Often, Triton X-100 seemed to work though the optimal reagent has to be selected for the respective protein and its hydrophilic|–phobic characteristics.

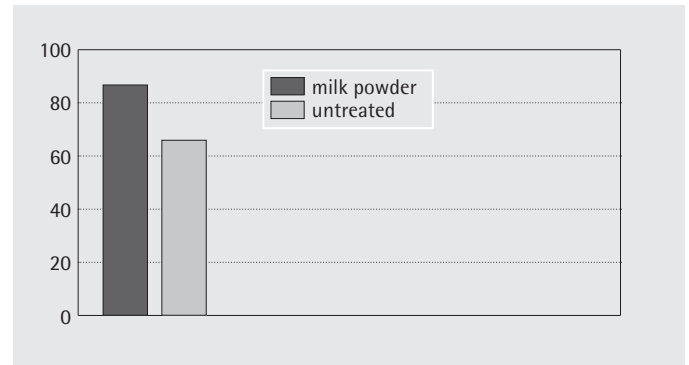

**Table 2:** Protein recovery (10 µg/ml BSA) with Vivaspin® PES 10 kDa after passivation

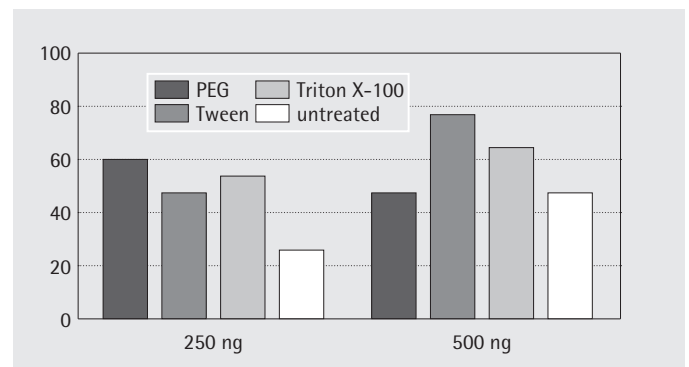

**Table 3:** Protein recovery (250 and 500 ng BSA) with Vivaspin® 2 PES 10 kDa after passivation

## Summary

Passivation is an appropriate method to achieve increasing sample recovery when using very dilute samples. In addition to skimmed milk, other proteins (BSA), detergents and compounds are possible. However, it should be noted that this is a general procedure, not specific for any particular application. Depending on the hydrophilic|–phobic character of the protein non-specific binding may be more or less of a problem and the suggested passivation solutions may lead to different results. Even with the Hydrosart membrane, which is recommended for dilute samples, passivation of the device will reduce losses on the plastic surface. One very important thing to remember is that the blocking agent is potentially introduced into the sample. It should be assured that this will not interfere with downstream analysis. For example, proteins must not be used for passivation if a pure protein is intended to be concentrated for x-ray crystallography, as even the smallest traces would interfere with the diffraction pattern. Other subsequent analyses methods include activity testing, gel electrophoresis or labelling are less problematic.

# Sales and Service Contacts

For further contacts, visit [www.sartorius-stedim.com](http://www.sartorius-stedim.com)

## Europe

### Germany

Sartorius Stedim Biotech GmbH  
August-Spindler-Strasse 11  
37079 Goettingen

Phone +49.551.308.0  
Fax +49.551.308.3289

[www.sartorius-stedim.com](http://www.sartorius-stedim.com)

Sartorius Stedim Systems GmbH  
Robert-Bosch-Strasse 5-7  
34302 Guxhagen

Phone +49.5665.407.0  
Fax +49.5665.407.2200

### France

Sartorius Stedim Biotech S.A.  
ZI Les Paluds  
Avenue de Jouques - BP 1051  
13781 Aubagne Cedex

Phone +33.442.845600  
Fax +33.442.845619

Sartorius Stedim France SAS  
ZI Les Paluds  
Avenue de Jouques - CS 71058  
13781 Aubagne Cedex

Phone +33.442.845600  
Fax +33.442.846545

### Austria

Sartorius Stedim Austria GmbH  
Franzosengraben 12  
1030 Vienna

Phone +43.1.7965763.18  
Fax +43.1.796576344

### Belgium

Sartorius Stedim Belgium N.V.  
Leuvensesteenweg, 248/B  
1800 Vilvoorde

Phone +32.2.756.06.80  
Fax +32.2.756.06.81

### Denmark

Sartorius Stedim Nordic A/S  
Hoerskaetten 6D, 1.  
2630 Taastrup

Phone +45.7023.4400  
Fax +45.4630.4030

### Hungary

Sartorius Stedim Hungária Kft  
Kagyló u. 5  
2092 Budakeszi

Phone +36.23.457.227  
Fax +36.23.457.147

### Italy

Sartorius Stedim Italy S.p.A.  
Via dell'Antella, 76/A  
50012 Antella-Bagno a Ripoli (FI)

Phone +39.055.63.40.41  
Fax +39.055.63.40.526

### Netherlands

Sartorius Stedim Netherlands B.V.  
Edisonbaan 24  
3439 MN Nieuwegein

Phone +31.30.6025080  
Fax +31.30.6025099

### Poland

Sartorius Stedim Poland Sp. z o.o.  
ul. Wrzesinska 70  
62-025 Kostrzyn

Phone +48.61.647.38.40  
Fax +48.61.879.25.04

### Russian Federation

000 „Sartorius ICR“ and 000 „Biohit“  
Uralskaya str. 4, Lit. B  
199155, Saint-Petersburg

Phone +7.812.327.5.327  
Fax +7.812.327.5.323

### Spain

Sartorius Stedim Spain SA  
C/Isabel Colbrand 10,  
Oficina 70  
Poligono Industrial de Fuencarral  
28050 Madrid

Phone +34.90.2110935  
Fax +34.91.3589623

### Switzerland

Sartorius Stedim Switzerland AG  
Ringstr. 24 a  
8317 Tagelswangen

Phone +41.52.354.36.36  
Fax +41.52.354.36.46

### U.K.

Sartorius Stedim UK Limited  
Longmead Business Park  
Blenheim Road, Epsom  
Surrey KT19 9 QQ

Phone +44.1372.737159  
Fax +44.1372.726171

## America

### USA

Sartorius Stedim North America Inc.  
5 Orville Drive, Suite 200  
Bohemia, NY 11716

Toll-Free +1.800.368.7178  
Fax +1.631.254.4253

### Argentina

Sartorius Argentina S.A.  
Int. A. Avalos 4251  
B1605ECS Munro  
Buenos Aires

Phone +54.11.4721.0505  
Fax +54.11.4762.2333

### Brazil

Sartorius do Brasil Ltda  
Av. Dom Pedro I, 241  
Bairro Vila Pires  
Santo André  
São Paulo  
Cep 09110-001

Phone +55.11.4451.6226  
Fax +55.11.4451.4369

### Mexico

Sartorius de México S.A. de C.V.  
Circuito Circunvalación Poniente No. 149  
Ciudad Satélite  
53100, Estado de México  
México

Phone +52.5555.62.1102  
Fax +52.5555.62.2942

## Asia | Pacific

### Australia

Sartorius Stedim Australia Pty. Ltd.  
Unit 5, 7-11 Rodeo Drive  
Dandenong South Vic 3175

Phone +61.3.8762.1800  
Fax +61.3.8762.1828

### China

Sartorius Stedim Biotech (Beijing)  
Co. Limited  
Airport Industrial Zone B  
No. 33 Yu'an Road  
Beijing 101300, Shunyi District

Phone +86.10.80426516  
Fax +86.10.80426580

Sartorius Stedim Biotech (Beijing)  
Co. Limited  
Shanghai Branch office  
Room 618, Tower 1, German Centre,  
Shanghai, PRC., 201203

Phone +86.21.28986393  
Fax +86.21.28986392.11

Sartorius Stedim Biotech (Beijing)  
Co. Limited  
Guangzhou representative office  
Room 704, Broadway Plaza,  
No. 233-234 Dong Feng West Road  
Guangzhou 510180

Phone +86.20.8351.7921  
Fax +86.20.8351.7931

### India

Sartorius Stedim India Pvt. Ltd.  
#69/2-69/3, Jakkasandra  
Kunigal Road, Nelamangala Tq  
Bangalore - 562 123

Phone +91.80.4350.5248  
Fax +91.80.4350.5253

### Japan

Sartorius Stedim Japan K.K.  
Kiba Park Bldg  
5-11-13 Kiba  
Koto-ku  
Tokyo 135-0042

Phone +81.3.5639.9981  
Fax +81.3.5639.9983

### Malaysia

Sartorius Stedim Malaysia Sdn. Bhd.  
Lot L3-E-3B, Enterprise 4  
Technology Park Malaysia  
Bukit Jalil  
57000 Kuala Lumpur, Malaysia

Phone +60.3.8996.0622  
Fax +60.3.8996.0755

### Singapore

Sartorius Stedim Singapore Pte. Ltd.  
1 Science Park Road,  
The Capricorn, #05-08A,  
Singapore Science Park 2  
Singapore 117528

Phone +65.6872.3966  
Fax +65.6778.2494

### South Korea

Sartorius Korea Biotech Co., Ltd.  
8th Floor, Solid Space B/D,  
PanGyoYeok-Ro 220, BunDang-Gu  
SeongNam-Si, GyeongGi-Do, 463-400

Phone +82.31.622.5700  
Fax +82.31.622.5799



◀▶ [www.sartorius-stedim.com](http://www.sartorius-stedim.com)